MagCore® Automated Nucleic Acid Extractor

MagCore® Automated Nucleic Acid Extractor Overview

		Standard Optional	Process Monitoring through your Smartphone MagCore® Plus //	Spectrophotometer Built-in MagCore® Super	NA Extraction and PCR Set up Mag Core® EDA
16 Sample	1-16 Samples			•	_
48 Sample	1-48 Samples				•
Spectrophotometer	Spectrophotometer	1// ' ' '		-	•
Touch Screen	Touch Screen		-	-	
UV	UV Decontamination		-	•	•
Barcode Sanner	Barcode Scanner		•	•	•
Thermo Phinter	Thermo Printer				
Upprakethvis (SS	Built-in Programs (Upgradeablevia USB ports, Plug&Play)		•		•
V	USB Output (USB flash drive not provided)		•		•
Mantaring	Progress Monitoring (Wireless)		•		
us	LIMS (Laboratory Information Management System)		•		
	HEPA Decontamination				
	Extraction+PCR set up				•
	PCRsetup				•


MagCore® Automated Nucleic Acid Extractors will keep you ahead





Automatic Optical Density Measurement

- Built-in spectrophotometer provides O.D. A₂₆₀ and A₂₈₀ measurement of individual samples. (O.D. detection range: ABS < 6.)
- A₃₂₀ Normalization
- Disposable cuvettes.

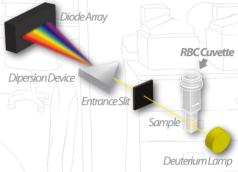

Thermo Printer and Barcode Scannel



Progress Monitoring





MagCore® Extractor System is a simple, fast and cost-effective instrument for automated purification of nucleic acids from a diverse range of sample sources. Featuring pre-programmed protocols and our unique magnetic-bead technology, MagCore System delivers efficient and consistent nucleic acid purification.

MagCore® Extractors are bench-top instruments ensuring efficient and cross-contamination free isolation of DNA/RNA. Built-in UV lamps allow to easily and efficiently decontaminate the instruments afterrun.

Flexibility

MagCore® Automated Extraction System allows you to save time without sacrificing consistency and purity. You can use one instrument to purify DNA and RNA from a broad variety of sample types: from blood to mouse tails and almost everything in between.

Ease Of Use

You will be provided with everything you need to run purifications, including pre-filled cartridges, specialized disposable tips and tubes. With the user-friendly interface and our user manuals, you are quaranteed to operate with ease.

Safety

MagCore® Automated Extraction System helps minimize cros-contamination by limiting hands-on procedures and tumaround time.

MagCore® System speeds the front-end processing, enabling you to do more tests in less time. And the Instrument is compact, soit can virtually fit into any lab.

Built-in Programs

All of our MaaCore Extractor models have built-in protocols for all of the kits we offer. Simply run the protocol by selecting the 3-digit codeprinted on the kit of interest.

Free upgrade of software and protocols can be downloaded from our website (www.rbcbioscience.com) and uploaded through the instrument RS232/USB ports.

Diverse Sample Purification

We offer extraction kits designed for Blood, Plasma, Cell, Tissue, FFPE Tissue and Plant samples, to fit all your research needs.

Competitive Price and Small Footprint

Easy To Use

Apply samples to instrument



Load Accessories



Select the number of the cartridge



Select Sample Volume



Push Start



A Beep Sound can be heard after protocol is completed












MagCore® Nucleic Acid Extraction Kits

MagCore® Automated Nucleic Acid Kits Selection Guide

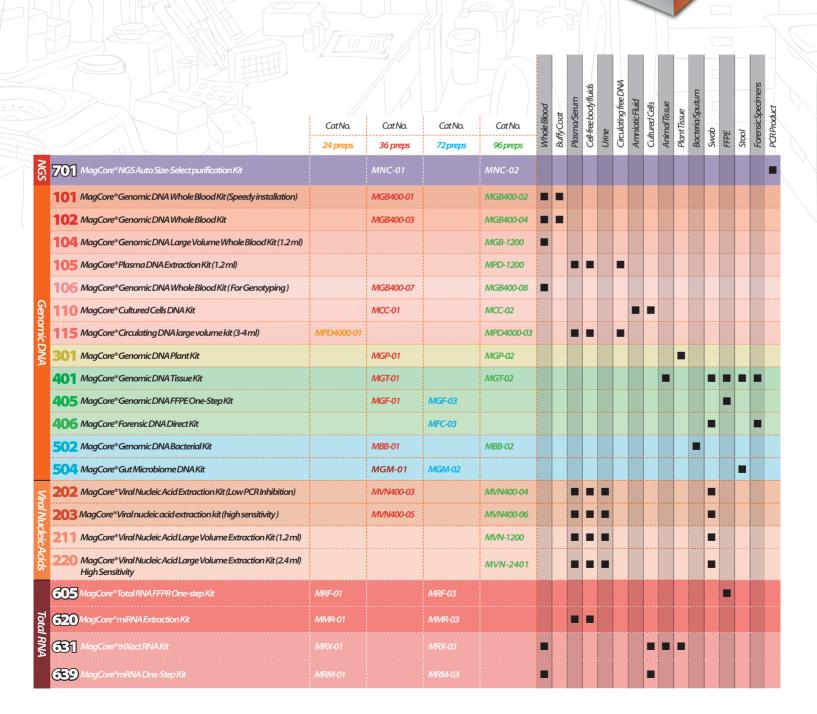


MagCore® Cartridge Design and Extraction Principle




MagCore® cartridges include all reagents needed for purifications, no additional handling is necessary. We minimize any possible contamination and spillage with an automated piercing step for our pre-sealed cartridges. RBC patented Heating Well and Separation Well in the cartridge provide a strong circular force to ensure efficient binding and washing during the extraction procedure.

MagCore® Nucleic Acid Extraction Kits contain all reagents and consumables needed for 36, 72 or 96 nucleic acid isolation reactions. The consumables consist of reagent cartridges, individually packaged tip sets, sample tubes and elution tubes.

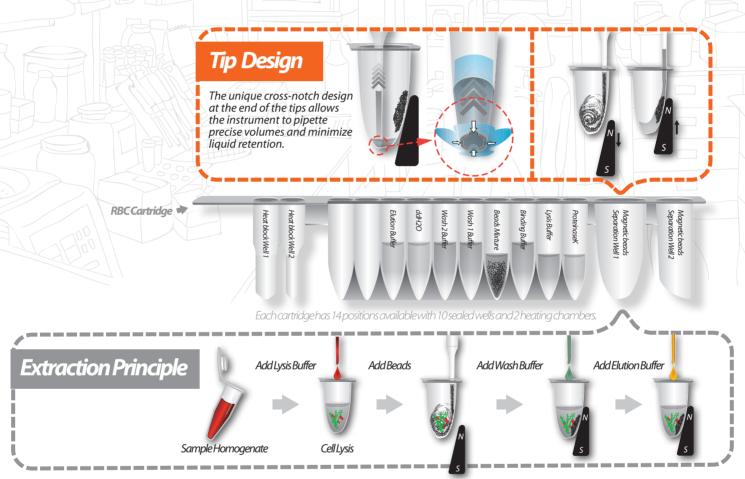

MagCore® Worldwide Patented Magnetic Beads

Our design: multiple core inside, cellulose coating.

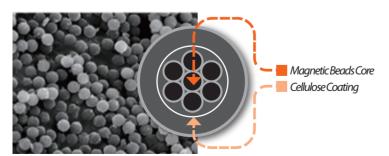

Particlerange	~20µm (inwater)
VSM	~40emu/g
Cellulosew%	~50%
Core size	~150nm
CoreVSM	~80emu/g
Core material	Fe ₃ O ₄
Bead Binding capacity	1 mg beads bind ~300µg Calfthymus DNA ~350µg human placenta DNA









MagCore® Automated Nucleic Acid Extractor

Full traceability and mobile monitoring on your smartphone

agCore® Plus ll



MagCore® Plus II is the newest robotic bench-top workstation for a fast and high-yield nucleic acid purification from virtually all molecular diagnostic biological, clinical and forensic sample types. With small footprint, light weight, user friendly interface, and a broad range of entirely built-in programs with free upgrades, 1-16 samples can be isolated simultaneously at your fingertip. The instrument simplifies your daily routine providina full traceability of kits and samples, through real-time mobile monitoring and a complete report that can be downloaded on a computer at the end of each run.



Worldwide Patented Magnetic Beads

Cellulose-coated magnetic beads, coupled with our patented binding and separation technology, quarantee high quality



Ideal for both DNA/RNA extraction

Built-in protocols are created for extracting nucleic acids from a wide range of samples, including whole blood, plasma (circulating free nucleic acid), tissue, bacteria, virus, plant and



Throughput up to 16 samples per run

From cartridge piercing to final eluate, all steps are performed by the instrument, that allows running 1 to 16 samples at one time, for a time-saving and flexible performance.



Full traceability of the samples and kits

A report in .csv format is generated at the end of each run and contains all relevant data: user's name, sample and kit barcode, protocol number, sample and elution volume, start and end time. The file, opened on a computer, can be subsequently processed by a LIMS.



Real-Time Mobile Monitorina

During the run, the instrument HMI can be accessed via Wi-Fi from your smartphone/tablet through our App, to see real-time information about the run processing status, remaining time and errors. Android and iOS compatible.



UVDecontamination

The equipped UV lamp minimizes the risk of crosscontamination and ensures user and product safety.



Built-in Programs (Upgradeablevia USB ports, Plug&Play)

MagCore® Plus II features built-in protocols for all the extraction kits we offer and is equipped with a USB port for free protocol and software upgrades.



Barcode Scanner

For sample and kit tracking and monitoring and an easier organization of the test results.



Easy To Use



Select User and Scan Barcodes



Select the code of the cartridae.



Select Sample Volume And Eluate Volume





A Beep Sound can be heard when the program completes



Open the run report on your computer







Samethroughput, smaller size




Barcode Scanner



Laboratory Information Management System (LIMS) Unidirectional LIMS device, Ethernet cable



Cartridge Design and Extraction Principle

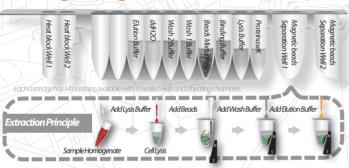

Mobile Monitoring with Android and iOS App



Specification

Model	Plus II	
System Method	Cellulose coated magnetic beads	
System Components	 Pipetting Unit: X and Y-axis movement for sample transfer and dispense. PLC module, HMI and Driver main board embedded in UV Light: power 8w, life duration 11,000 hrs Heating Block: RT-90°C Display Screen: 7-inch color touch panel Accessories: T-racks, cartridge racks, barcode scanner, waste box 	
Power Supply	Voltage: AC 100V~240V; Frequency: 50/60Hz	
Dimension	W523 x D602 x H605 (mm) / W21 x D23.7 x H23.8 (inches)	
Net Weight	70kg/154.35lbs	

Operating Parameters

Operating Environment

3 , ,	1-16 samples per batch	Temperatures allowed during transportation, storage, and packaging	
Processing Time	30-90 minutes (depends on sample type and method)		
	200 µl/400 µl/1,200 µl/ 4ml * depending on the program.		
Elution Volume	30µl/40µl/60µl/100µl/150µl/200 µl * depending on the program.	Temperatures allowed during operation	18℃-30°C
Yield	Average 6µg Genomic DNA from 200µl human whole blood	Pollution Degree	Level 2
Purity	DNA: O.D A ₂₆₀ / ₂₈₀ ratio 1.8 ± 0.1 RNA: O.D A ₂₆₀ / ₂₈₀ ratio 2.0 ± 0.2		
	30-60µl 20%; 60-100µl 10%; 100-1000 µl 4%		













Automatic Extraction and Smart PCR Setup



Worldwide Patented Magnetic Beads

Cellulose-coated magnetic beads, coupled with our patented binding and separation technology, quarantee high quality extracts.

Ideal for both DNA/RNA extraction

Built-in protocols are created for extracting nucleic acids from a wide range of samples, including whole blood, plasma (circulating free nucleic acid), tissue, bacteria, virus, plant and forensic.

Automatic Optical Measurements of OD Values

Built-in spectrophotometer and our optical module provide users the option to automatically measure OD values and concentration of final eluates upon completion of the nucleic acid extraction process.

WDecontamination

The equipped W lamp minimizes the risk of cross-contamination and ensures user and product safety.

HEPA Decontamination

HEPA system is installed in the PCR set up zone to eliminate contamination and ensure the purity of the samples.

Automatic Barcode Scanning

Barcode reading of primary tube simplifies the tracking of the samples and ensures traceability.

Throughput up to 48 samples

From cartridge piercing to final elute, all steps are performed by the instrument, that allows running 1-48 samples. The instrument processed samples in batches of 24 at a time and has run setting for 24/48 sample to be processed at once, for time saving and flexible performance.

Smart PCR setup

After pre-mix buffer configuration is set, the instrument automatically calculates the required volumes of the primers and perform liquid level detection. Pre-run report is generated to avoid errors and énsure optimal performance.







Automatic Optical Density Measurement

- Built-in spectrophotometer provides O.D. A230, A₂₆₀ and A₂₈₀ measurement of individual
- (O.D. detection range: ABS < 6.)
- A₃₂₀ Normalization
- Disposable cuvettes.



EDA System Connectivity

With LIMS compatibility EDA system can be connected to various thermocyclers and laboratory computer to set up automatic information exchange and analysis workflow.



SmartPCRSetUp

Zone B provides automatic PCR set up which is compatible with different thermocylers for the further process. The system has 10 SBS positions, option of keeping premixes cool and the option of using 0.1 and 0.2 ml standard tubes, along from customised tubes from other PCR supplier.



'	
Model	EDA /
System Method	Cellulose coated magnetic beads
System Components	1. Pipetting Unit: dispensing, transferring, 2 X-Y axis model 2. Electric Control: PLC model and Arm-based main board embedded in 3. UV Light: power 8W, life duration 11,000hrs 4. Contamination Prevention: HEPA filtration, Pipette tips with filter 5. Heating Block: RT-959C 6. Nucleic Acid Concentration Detection Source: D2 lamp 7. Wavelength Detection: 230nm, 260nm, 280nm 8. Liquid detection sensor: PCR setup Zone 9. Automatic Barcode Scanning: Primary tube 10. Thermoelectric Cooler 4~10°C 11. Accessories: T-rack, Cartridge racks, trash drawer, 96-well PCR plate rack, 32-well 1.5ml tube rack 12. Stand (optional)
Power Supply	Voltage: AC 200-240V; Frequency: 50/60Hz
Dimension	W1240*D920*H830(mm)/W48.81 x D36.22 x H32.67(inches)
Net Weight	335 Kg / 739 lbs
Net Weight (with stand)	460 Kg / 1,014 lbs

Operating Parameters

Processing Capacity 1-48 samples (in batches of 24)	
Processing Time	30-120minutes (depends on sample type)
Sample Volume	200/400/1200/4000μl (application dependent)
Elution Volume	30/40/60/100/150/200 μl (application dependent)
Yield	200ul whole blood (average 6ug gDNA) 400ul whole blood (average 12ug gDNA)
Purity	DNA: O.D A ₂₆₀ /A ₂₆₀ ratio 1.8±0.1 RNA: O.D A ₂₆₀ /A ₂₆₀ ratio 2.0±0.2
Pipetting Accuracy	Extraction Zone: 40µl <5%; 60µl <2%; 100-900µl <1.5% PCR setup Zone:

Operating Environment

i

Ī

Processing Capacity	1-48 samples (in batches of 24)		
Processing Time	30-120minutes (depends on sample type)	Temperatures allowed during transportation, storage, and packaging	15°C-35°C
Sample Volume	200/400/1200/4000 µl (application dependent)	transportation, storage, and packaging	
Elution Volume	30/40/60/100/150/200 µl (application dependent)	Temperatures allowed during operation	18℃-30°C
Yield	200ul whole blood (average 6ug gDNA) 400ul whole blood (average 12ug gDNA)	Pollution Degree	Level 2
Purity	DNA:O.D A ₂₆₀ /A ₂₆₀ ratio 1.8±0.1 RNA:O.D A ₂₆₀ /A ₂₆₀ ratio 2.0±0.2		
Pipetting Accuracy	Extraction Zone: 40µl <5%; 60µl <2%; 100-900µl <1.5% PCR setup Zone: CV<1%, 5~50µl; CV<5%, 2~5µl; CV<10%, 1µl		











The first instrument with a built-in Spectrophotometer

MagCore® Super















Automatic Optical Measurements of OD Values

Built-in spectrophotometer and our optical module provide users the option to automatically measure OD values and concentration of final eluates upon completion of the nucleic acid extraction process.

Test Report

Test results can be saved in the instrument, downloaded through he USB port and/or printed by the thermal printer. USB Output (USB flash drive not provided)

USB Output allows users to conveniently save tests reports in excel format and upload system updates with a USB flash drive.

Thermal Printer

Test reports are available in hard copy.

Laboratory Information Management System (LIMS)

Test results are automatically saved after optical measurements. You can save up to 1,600 tests in LIMS. Data can be easily transferred to a printer or computer in the same network and the report file can be subsequently processed by a LIMS.



Worldwide Patented Magnetic Beads

Cellulose-coated magnetic beads, coupled with our patented binding and separation technology, quarantee high quality extracts.



Ideal for both DNA/RNA Extraction

Built-in protocols are created for extracting nucleic acids from whole blood, plasma, tissue cell, plant cell, bacteria cell and virus samples.



UVDecontamination

The equipped UV lamp minimizes the risk of cross-contamination and ensures user and product safety.



High Capacity of 16 Samples

The instrument and protocols allow running up to 16 samples at one time, providing time-saving and flexible operation.



Built-in Programs (Upgradeable via USB ports, Plug&Play)

MagCore Super has built-in protocols for all of the extraction kits we offer. Simply run the protocol by selecting the 3-digit code printed on the kit of interest. MagCore® is equipped with a USB port. Free upgrade of software or protocols can be downloaded from our website (www.rbcbioscience.com).



Touch Screen with User-Friendly Interface

An integrated 7-inch full-color touch screen with user-friendly interface offersease in operation. Only one touch is required to run your daily work.



Barcode Scanner

It enables sample tracking and monitoring throughout the entire purification process and helps organize test results.



Progress Monitoring

Remote (wireless) HMI device automatically transfers the data to you Android smartphone.



Load Samples



Install Accessories



Select the code of the cartridge.



Select Sample Volume



Press Start




A Beep Sound can be heard after the program is complete!








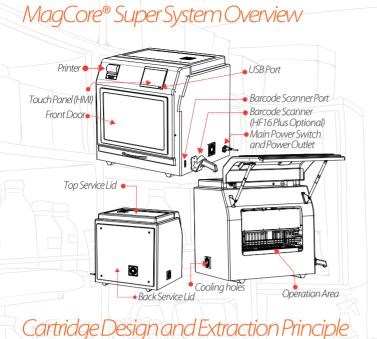

Automatic Optical Scanning and Retrieval of ODValues

 The optical module provides O.D. A₂₆₀ and A₂₆₀ measurement of individual samples.(O.D. detection range: ABS < 6.)

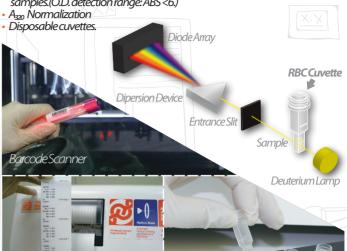





Process Monitoring (Wireless-Android Only) (optional)



Specification

Add Lysis Buffer

Add Beads

Add Wash Buffer Add Elution Buffer

Heat block Well

Model	Super /		
System Method			
System Components	 Pipetting Unit: X and Y-axis movement for sample transfer and dispense. Electric Control: PLC module and ARM-based main board embedded in UV Light: power 8w, life duration 11,000hrs Heating Block: RT-90° C OD Detection Rouge: ABS 0 - 2.5 Detection Source: D2 lamp Detection Wavelength: 260nm, 280nm Display Screen: 7-inch color touch panel Accessories: T-racks, cartridge racks, cuvette racks, barcode scanner, thermal printer 		
Power Supply	Voltage: AC 100V~ 240V; Frequency: 50/60Hz		
Dimension	W760xD700xH770(mm)/W29.92xD27.55xH30.31 (inches)		
Net Weight	78kg/171.99lbs		

Operating Parameters

_	
Processing Capacity	1-16 samples per batch
Processing Time	30-90 minutes (depends on ample type and method)
Sample Volume	200 µl/400 µl/1,200 µl /4ml * depending on the program.
Elution Volume	30µl/60µl/100µl/150µl/200µl
Yield	Average 6µg Genomic DNA from 200µl human whole blood
Purity	DNA: O.D $A_{260}/_{280}$ ratio 1.8 ± 0.1 RNA: O.D $A_{260}/_{280}$ ratio 2.0 ± 0.2
Pipetting Accuracy	500µl ≤ 4%

Operating Environment

Temperatures allowed during transportation, storage, and packaging	15°C-35°C
 Temperatures allowed during operation	18℃-30°C
Pollution Degree	Level 2













For selective DNA purification and size selection for NGS

MagCore® NGS Auto Size-Select purification Kit is designed for fast DNA fragment cleanup and size selection. The cleanup program results in 80% or higher recovery while removing all adapter dimers in sample that is then applicable during the preparation of Next Generation Sequencing (NGS) library. For DNA fragment size selection, the program selectively purifies nucleic acid fragments from 100-200 bp or from 200-500 bp in just minutes which is also suitable for library preparation. The entire purification procedure is fully automated and can be performed in 35 minutes.

701

MagCore®NGS Auto Size-Select purification Kit

Magnetic beads
Separation Well 2
Separation Well 1
Separation Well 1
Separation Well 1
Separation Well 1
Seads Mixture 1000ul
Wash Buffer 1 1000ul
Wash Buffer 1 1000ul
Wash Buffer 1000ul
Button Buffer 1000ul
Button Buffer 1000ul
Heat block Well 1
Heat block Well 1

Features

- 1. Automated DNA Cleanup and Size Selection ~ 30min
- 2. Precise Size Range Selection
- 3. High Recovery Rate
- 4. Easy and Rapid Workflow
- 5. Consistent and Reliable

Applications

The MagCore® NGS Auto Size-Select purification Kit is designed to purify DNA for various downstream applications, including:

- 1. DNA library peparation for NGS
- 2. PCRorreal-time PCR




Product Introduction and Average Recovery

Program	Program Description	Avarage Recovery (%)
701A	DNA Cleanup	>80%
701B	DNA Size Selection 200-500bp	40~60%
701C	DNA Size Selection 100-200bp	>50%

To meet the requirement of NGS, it is better to produce a NGS library with DNA fragment sizes between 100~200 bp(https://doi.org/10.1016/B978-0-12-418687-3.00015-X) or 200~500 bp (https://doi.org/10.1007/s11262-018-1568-x).


Comparison of DNA Cleanup methods

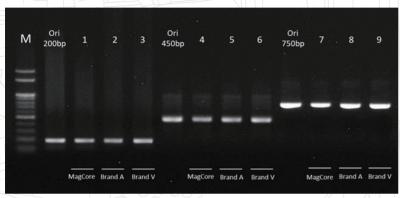

M: 100bp Marker. Lane 1, 4, 7: MagCore NGS Auto Size-Select purification Kit cleanup DNA sample with 701A program. Lane 2, 5, 8: Brand A manually cleanup DNA sample. Lane 3, 6, 9: Brand V manually cleanup DNA sample.

MagCore NGS Auto Size-Select purification Kit: results of the automatic cleanup of DNA sample are comparable with Brand A and Brand V manual kits, resulting in high purity product with >80% recovery.

Human genomic library prepared with DNA Library Prep Kit was analyzed by an Agilent Technologies 2100 Bioanalyzer

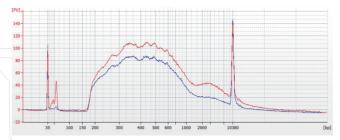

Red Line: Before DNA Clean-up Blue Line: After DNA Clean-up

Precise Size Selection



M:100bpMarker. Lane 1:MagCore NGS Auto Size-Select purification Kit sizeselect 200-500bp with 701B program. Lane 2:Brand A manually size select 200-500bp. Lane 3:Brand V manually size select 200-500bp.

MagCore NGS Auto Size-Select purification Kit is completely aotomated and can effectively size select DNA from 200-500bp, while Brand A and Brand V contain larger DNA fragments in the final product.

MagCore®Automated Nucleic Acid Kits Specification

MagCore®Super/HF16Plus/PlusII/EDA

Cat No.	CatNo.	Di vanina Timo
36 preps	96preps	Running Time

Wageore Til 107 Compact							
	Cat No.	Di maina Tima					
	96preps	RunningTime					

MaaCoro® HE16 /Compact

701

Cartridge Code

MagCore®NGS Auto Size-Select purification Kit

For 50 µl sample volumes.

Contents: Pre-Filled Cartridges, Disposable Tip & Holder Sets, Sample Tubes, Elution Tubes

Shelf life: 18 months

İ						
	MNC-01	MNC-02	35 min	MNC-01	MNC-02	35 min

Cat No.

36 preps









MagCore® Nucleic Acid Extraction Kits



For extraction of genomic DNA from human whole blood sample

MagCore® Genomic DNA Whole Blood Kit is designed for extraction of total DNA (including genomic, mitochondrial and viral DNA) from whole blood, plasma, serum and buffy coat. The pre-filled cartridge contains proteinase K, a chaotropic salt, and guanidine hydrochloride for cell lysis and protein degradation. They enhance the binding between cellulose-coated magnetic beads and DNA. High quality DNA is eluted by low salt elution buffer or water after the removal of contaminants. Purified DNA of approximately 20-30 kb is suitable for PCR or other downstream applications.


101

MagCore® Genomic DNA Whole Blood Kit (Speedy installation)



102

MagCore® Genomic DNA Whole Blood Kit

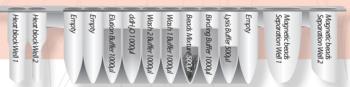

104

MagCore® Genomic DNA Large Volume Whole Blood Kit (1.2 ml)



106

MagCore® Genomic DNA Whole Blood Kit (For Genotyping)



l-eatures

- 1. High performance of purified DNA in downstream applications such as gPCR.
- 2. High analytical sensitivity
- 3. Cartridges are pre-filled and sealed to prevent contamination.
- 4. Walkaway processing improves work efficiency.

Applications

The MagCore® Genomic DNA Whole Blood Kit is designed to allow automated processing of multiple sample types in the same run. Sample types include:

- 1. Fresh and frozen whole blood
- 2. Buffycoat
- 3. Bodyfluids


High quality DNA available for various downstream applications, including:

- 1. PCR and real time PCR
- 2. Genotyping or sequencing
- 3. SNP,STR




DNA quality unaffected by anticoagulants

Anticoagulant	A ₂₆₀ / ₂₈₀	Conc.(ng/µl)
1.EDTA	1.87	99.3
2. Sodium Citrate	1.87	86.1
3. Lithium Heparin(4ml)	1.95	106.02
4. Sodium Fluoride	1.91	113.16
5. Lithium Heparin(9ml)	1.95	101.34

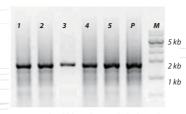


Consistent yield and purity

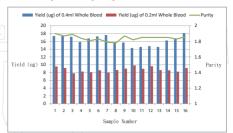

It is shown that there is a scalability of DNA isolation from blood samples (20, 50, 80, 100, 150, and 200 µl) by the MagCore® HF16 The amount of DNA was determined by A_{200} measurement.





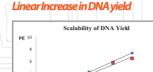

MagCore® Automated Nucleic Acid Kits Specification

MagCore®Super/HF16Plus/PlusII/EDA

MagCore® HF16 / Compact / HF48

	_\C
Cartridge Code	9-1-
	36

36 preps	96preps
CatNo.	Cat No.

Running Time

CatNo. CatNo. 36 preps 96preps

Running Time

101

MagCore®Genomic DNA Whole Blood Kit (Speedy installation)

For 200 and 400 µl sample volumes.

Contents: Pre-Filled Cartridges (Including Proteinase K), Disposable Tip & Holder Sets, Sample Tubes, Elution Tubes

Shelf life: 12 months

MGB400-01

39min (samplevolume:200 µl) 50min (samplevolume:400 µl)

MGB400-01

MGB400-02 44 n

44 min (sample volume:200 μl) 57 min (sample volume:400 μl)

102

MagCore®Genomic DNA Whole Blood Kit

MGR400-02

MGB400-04

For 200 and 400 µl sample volumes.

Contents: Pre-Filled Cartridges (Including Proteinase K), Disposable Tip & Holder Sets, Sample Tubes, Elution Tubes

Shelf life: 18 months

MGB400-03

39min (sample volume:200 µl) 50min (sample volume:400 µl)

MGB400-03

MGB400-04

44 min (sample volume:200 µl) 55 min (sample volume:400 µl)

104

MagCore®Genomic DNA Large Volume Whole Blood Kit (1.2 ml)

For 1200 µl sample volumes.

Contents: Pre-Filled Cartridges (Including Proteinase K), Disposable Tip & Holder Sets, Sample Tubes, Elution Tubes

Shelf life: 18 months

MGB-1200

76min (samplevolume:1200 µl)

MGB-1200

83 min (sample volume:1200 µl)

106

MagCore® Genomic DNA Whole Blood Kit (For Genotyping)

For 200 and 400 µl sample volumes.

Contents: Pre-Filled Cartridges (Including Proteinase K), Disposable Tip & Holder Sets, Sample Tubes, Elution Tubes

Shelf life: 18 months

MGB400-07

MGB-400-08

41 min (samplevolume:200 μl) 53 min (samplevolume:400 μl) * optical detection is not provided

MGB400-07

MGB-400-08

44 min(samplevolume:200 μl) 57 min(samplevolume:400 μl) (Applicable models: HF16, Compact)

Enzyme Selection Guide

Product	Contents	Cat.No.
Proteinase K Set	11 mg Proteinase K,1.25 ml PK Storage Buffer	PK011
04/ 4	50µl RNaseA (50mg/ml)	RN050
KNASE A	130 µIRNase A (50mg/ml)	RN130









RBC Bioscience Corp.
www.rbcbioscience.com
info@rbcbioscience.com





For extraction of circulating free DNA from human plasma or serum

MagCore® Plasma DNA Extraction Kit is designed for purifying circulating free DNA from human serum or plasma using MagCore® automated extraction systems. The kit contains all required reagents and labware for automated purification based on magnetic-particle technology. Reagents necessary for a complete process are supplied and pre-filled in the cartridges, which can be easily loaded into the MagCore®instrument.

105

MagCore® Plasma DNA Extraction Kit (1.2 ml)

Magnetic beads
Separation Well 1

Separation Well 1

Empty
Lysis Buffer 1200µl

Binding Buffer 1500µl

Beads Mixture 500µl

Mash 1 Buffer 1000µl

Mash 2 Buffer 1000µl

DEPC Water 1000µl

DEPC Water 1000µl

DEPC Water 1000µl

Empty

Empty

Empty

Heatblock Well 1

Features

- 1. High performance of purified DNA in downstream applications such as gPCR.
- 2. Efficient recovery of fragmented DNA
- 3. Pre-filled and sealed buffer cartridges prevent contamination.
- 4. Nophenolorchloroform extraction
- 5. Efficient removal of contaminants and inhibitors.

Applications

The MagCore® Plasma DNA Extraction Kit efficiently purifies circulating free DNA. Sample types include:

- 1. Human plasma
- 2. Human serum

High quality DNA available for various downstream applications, including:


- 1. PCR and real-time PCR
- 2. Next Generation Sequencing (NGS)



High-quality of cell-free DNA from plasma and serum

Quality and quantity analysis of the cell-free DNA by Agilent Bioanalyzer 2100. Superior quality DNA are available from plasma and serum samples by using MagCore® products.


Sample ID	Sample type	Target Conc. (pg/μl)	35 bp marker Conc. (pg/µl)	10380 bp marker Conc. (рд/µl)
1	Human plasma	77.94	125	<i>7</i> 5
2	Human serum	138.48	125	<i>7</i> 5

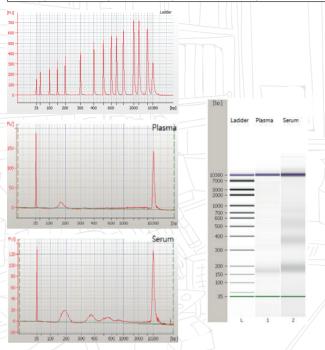

High stability of cell-free DNA

Quality and quantity analysis of the cell-free DNA by Agilent 2100 Bioanalyzer.It is shown by the evidence that cell-free DNA of high quality and stability is possible to extract from 5 different plasma samples by using MagCore® procudts.

Sample ID	Sample type	Target Conc.(pg/μl)	35 bp marker onc. (pg/μl)	10380 bp marker Conc. (рд/µl)
1	Human plasma	65.95	125	<i>7</i> 5
2	Human plasma	<i>7</i> 6.63	125	<i>7</i> 5
3	Human plasma	87.22	125	<i>7</i> 5
4	Human plasma	57.96	125	<i>7</i> 5
5	Human plasma	60.53	125	<i>7</i> 5

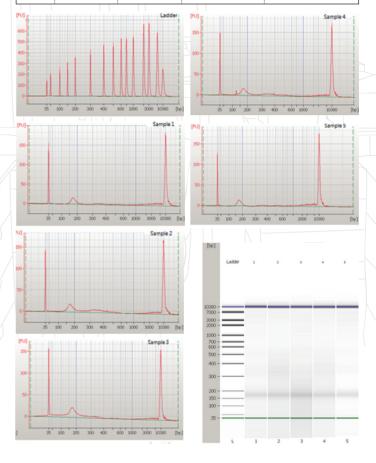

MagCore® Automated Nucleic Acid Kits Specification

MagCore®Super/HF16Plus/PlusII/EDA			MagC	ore® HF16/C	Compact/HF48
Cartridge Code	CatNo.	Running Time		CatNo.	Running Time
Carmage Code	96preps			96preps	Kuririing III'le

MagCore® Plasma DNA Extraction Kit (1.2 ml)

For 1200µl sample volumes. Contents: Pre-Filled Cartridges, Proteinase K, PK Storage Buffer, Disposable Tip & Holder Sets, Sample Tubes, Elution Tubes Shelf life: 18 months

|--|

Enzyme Selection Guide

Product	Contents	Cat.No.
Proteinase K Set	11 mg Proteinase K, 1.25 ml PK Storage Buffer	PK011
RNase A	50 µl RNase A (50mg/ml)	RN050
	130μlRNaseA (50mg/ml)	RN130













For genomic DNA extraction from cultured cells and amniotic fluid

MagCore® Cultured Cells DNA Kit is designed to extract genomic DNA from up to 5x106 cultured cells using MagCore® automated extraction systems. The kit contains all required reagents and labware for automated extraction using magnetic-particle technology. Reagents are supplied in prefilled cartridges, which can be easily loaded into the MagCore® instrument.

MagCore® Cultured cells DNA Kit

Heat block Well :


Features

- 1. High performance of purified DNA in downstream applications such as gPCR.
- 2. Efficient isolation of DNA from up to 5 x 106 cells.
- Cartridges are pre-filled and sealed to prevent contamination.
 No phenol or chloroform extraction.
- 5. Efficient removal of contaminants and inhibitors.

Applications

High quality DNA available for various downstream applications, including:

- 1. PCR and real-time PCR
- 2. Next Generation Sequencing (NGS)

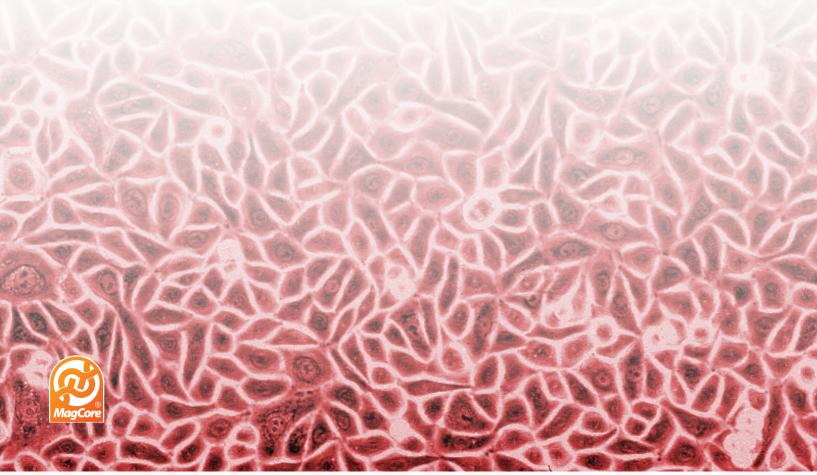


DNA quality analysis by Spectrophotometer

DNA of three different cultured cells were extracted by the employment of MagCore® HF16 and MagCore® Cultured Cells DNA Kit (200µl sample volume). The following data show the yield and purity of DNA of three samples.

Sample 1:COS7 (1.2x10⁷) Sample 2: Junkat (1x10°) Sample 3: HEK293 (1x10°)

	1	2	3
Yield (µg)	16.44	13.08	24.42
DNA purity (A ₂₆₀ / ₂₈₀)	1.93	1.87	1.92

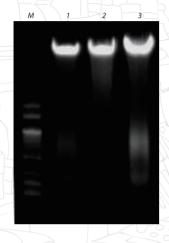

Amniotic fluid DNA extration

DNA quality is checked by the NanoDrop spectrophotometer and agarose gel electrophoresis after DNA purification from amniotic-fluid cells. Cells were harvested from 10~15 ml amniotic fluid samples at pregnancy weeks 16-18 by centrifugation for 10 minutes at 3000 rpm.

SampleID	Conc. (pg/µl)	A ₂₆₀ / ₂₈₀	A ₂₆₀ / ₂₃₀	Yield (μg)
(1) GPT2967P	22.68	1.84	1.07	1.36
(2) GPT2952P	33.98	1.92	1.57	2.04
(3) NP679P	79P 33.65		1.34	2.02
(4) NP777P	NP777P 39.22 1.86		1.20	2.35
(5) MP795P	22.05	1.79	0.97	1.32

by NanoDrop ND-1000

Table 1. DNA quality analysis by NanoDrop.

Sample ID (1) to (5) are different amniotic fluid samples. It is shown that MagCore® HF16 system can purify $1\sim2\,\mu g$ DNA from $10\sim15\,m$ amniotic fluid samples and DNA purity is with the $A_{200}/_{200}$ ratio of around 1.8±0.1



Figure 1. Genomic DNA isolations were run by gel electrophoresis on 1% agarose gel.

Lane M: RBC 1kb ladder marker

Lane 1:GPT2967P

Lane 2: GPT 2952P

Lane 3: NP679P Lane 4: NP777P

Lane 5: NP795P

MagCore® Automated Nucleic Acid Kits Specification

MagCore®Super/HF16Plus/PlusII/EDA

MagCore® HF16 / Compact / HF48

		. 3			- 5 -		, , , , , , , , , , , , , , , , , , ,
Cartridge Code	CatNo.	CatNo.	Running Time	Cat No.		CatNo.	Running Time
	36 preps	96preps		36 preps		96preps	

MagCore®Cultured Cells DNA Kit

For 200µl sample volumes (up to 5x10⁶ cells)
Contents: Pre-Filled Cartridges, Proteinase K, PK Storage Buffer, Disposable Tip & Holder Sets, Sample Tubes, Elution Tubes Shelf life: 18 months

 $39 \min (sample volume: 200 \mu l, up to <math>5x \cdot 10^6 cells)$ MCC-01SP MCC-02SP MCC-01 MCC-02 44 min (sample volume: 200 μl, up to 5 x 10 cells)

Enzyme Selection Guide

Product	Contents	Cat.No.
Proteinase K Set	11 mg Proteinase K, 1.25 ml PK Storage Buffer	PK011
RNase A	50 µl RNaseA (50mg/ml)	RN050
	130 µl RNase A (50mg/ml)	RN130











For extraction of cf DNA from 4ml plasma or serum

MagCore® Circulating DNA large volume kit is designed for purifying circulating free DNA from human serum or plasma using MagCore® automated extraction systems. The kit contains all required reagents and labware for automated purification based on magnetic-particle technology. Reagents necessary for a complete process are supplied and pre-filled in the cartridges, which can be easily loaded into the MagCore®instrument.

115

MagCore® Circulating DNA large volume kit (4ml)

Magnetic beads
Separation Well 2

Magnetic beads
Separation Well 1

Binding Buffer 1800µl

Lysis Buffer 1800µl

Beads Mixture 500µl

Beads Mixture 1000µl

Mash 1 Buffer 1000µl

Mash 2 Buffer 1000µl

AdH,O 1000µl

ddH,O 1000µl

ddH,O 1000µl

Lysis Buffer 1000µl

ddH,O 1000µl

Heatblock Well 1

Features

- 1. High performance of purified DNA in downstream applications such as qPCR.
- 2. Efficient recovery of fragmented DNA
- 3. Pre-filled and séaled buffer cartridges prevent contamination.
- 4. No phenol or chloroform extraction
- 5. Efficient removal of contaminants and inhibitors.

Applications

The MagCore® Plasma DNA Extraction Kit efficiently purifies circulating free DNA. Sample types include:

- 1. Human plasma
- 2. Human serum

High quality DNA available for various downstream applications, including:

- 1. PCR and real-time PCR
- 2. Next Generation Sequencing (NGS)



High stability of cell-free DNA

Quality and quantity analysis of the cell-free DNA by Agilent 2100 Bioanalyzer.It is shown by the evidence that cell-free DNA of high quality and stability is possible to extract from 5 different healthy people plasma samples by using MagCore products.

Sample ID	Sample type	Plasma (ml)	Qubit (ng/μl)	Ctvalue (GAPDH gene)	Bioanalyzer Conc. (pg/µl)	35 bp marker Conc. (pg/µl)	10380 bp marker Conc. (pg/µl)	
1	Human plasma	4	0.262	18.52	127.21	125	75	
2	Human plasma	4	0.250	18.59	124.62	125	75	
3	Human plasma	4	0.220	18.78	111.11	125	<i>7</i> 5	
4	Human plasma	4	0.290	18.45	153.24	125	75	
5	Human plasma	4	0.278	18.49	130.97	125	<i>75</i>	


Sample 1 Sample 2 Sample 3 Sample 4 Sample 5

Comparison of cell-free DNA from 1.2 ml and 4 ml plasma

Quality and quantity analysis of the cell-free DNA by Agilent Bioanalyzer 2100. Superior quality DNA are available from plasma samples by using MagCore products.

Sample ID	Kits Code	Sample type	Plasma (ml)	Qubit (ng/μl)	Ctvalue (GAPDH gene)	Bioanalyzer (pg/μl)	35 bp marker Conc. (pg/μl)	10380 bp marker Conc. (pg/µl)
1	105	Human plasma	1.2	0.061	23.86	24.70	125	<i>7</i> 5
2	105	Human plasma	1.2	0.058	23.96	22.98	125	<i>7</i> 5
3	115	Human plasma	4	0.203	20.04	117.0	125	<i>75</i>
4	115	Human plasma	4	0.194	20.26	104.7	125	75

Code 105:MagCore® Plasma DNA Extraction Kit Code 115:MagCore® Circulating DNA large volume kit

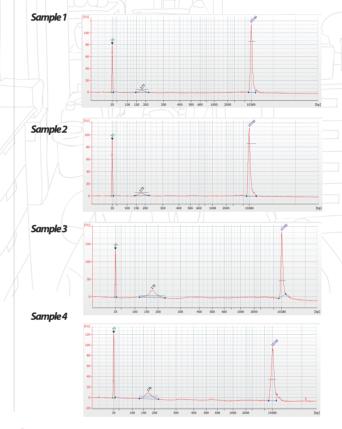

MagCore® Automated Nucleic Acid Kits Specification

MagCore®Super/HF16Plus/PlusII/EDA					MagCore® HF16 / Compact			
CartridaoCodo	CatNo.	Cat No.	Running Time	Cat No.		CatNo.	Running Time	
Cartridge Code	24 preps	96preps		24 preps		96preps	Kunning iime	

MagCore® Circulating DNA large volume kit

For 4000µl sample volumes

Contents: Pre-Filled Cartridges, Proteinase K, PK Storage Buffer, Disposable Tip & Holder Sets, Sample Tubes, Elution Tubes Shelf life: 18 months

MPD4000-01		MPD4000-03	146 min (sample volume:4000 µl) *ontical detection is not provided	MPD4000-01	MPD4000-03	146min (sample volume:4000 µl)
------------	--	------------	--	------------	------------	--------------------------------

Enzyme Selection Guide

Product Contents		Cat.No.
Proteinase K Set	11 mg Proteinase K, 1.25 ml PK Storage Buffer	PK011
	50µl RNaseA (50mg/ml)	RN050
RNase A	130 µI RNase A (50mg/ml)	RN130









RBC Bioscience Corp. www.rbcbioscience.com info@rbcbioscience.com





For extraction of viral DNA/RNA from serum, plasma and cell-free body fluids

The MagCore® Viral Nucleic Acid Extraction Kit is designed for purifying viral DNA and RNA from serum, plasma, and cell free body fluids. MagCore® magnetic particle technology delivers high-quality DNA/RNA that is suitable for direct use in downstream applications such as amplification or other enzymatic reactions. To minimize the risk of cross-contamination, plastic consumables inside the Kit are DNase/RNase treated and the operation system is designed to individually process samples at the same time. Multiple protocols are installed on the instrument and optimized for different sample volumes.

202

MagCore® Viral Nucleic Acid Extraction Kit (Low PCR Inhibition)



211

MagCore® Viral Nucleic Acid Large Volume Extraction Kit (1.2 ml)



Features

- 1. High performance of purified DNA in downstream applications such as qPCR.
- 2. High analytical sensitivity.
- 3. Cărtridges are pre-filled and sealed to prevent contamination.
- 4. Walkaway processing improves work efficiency.

Applications


MagCore® Viral Nucleic Acid Extraction Kit is designed to allow automated processing of multiple sample types in the same run. Sample types include:

- 1. Cell-free body fluids
- 2. Plasma and serum
- 3. CSF
- 4. Urine



HBV detection by Real-Time PCR

The isolations of HBV nucleic acid from samples containing different concentrations of HBV were subsequently detected by Real-Time PCR

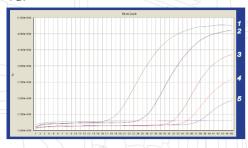

Sensitivity test for HBV detection

Lane 1	1x10°HBV serums.
Lane 2	1x10 ¹ HBV serums.
Lane 3	1x10 ² HBV serums.
Lane 4	1x10³HBV serums.
Lane N	Negative control.
Lane M	RBC 100 bp DNA Ladder Marker.

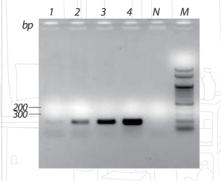


Figure 1. Nested PCR results of HBV at different concentrations. Viral nucleic acids were purified from samples containing different amount of HBV using MagCore® Viral Nucleic Acid Extraction Kit.

Sensitivity test for HCV detection

Lane 1	5x10 ^o HCV serums.
Lane 2	5x10 ¹ HCV serums.
Lane 3	5x10 ² HCV serums.
Lane 4	5x10³HCV serums.
Lane 5	5x10⁴HCV serums.
Lane 6	5x10⁵HCV serums.
Lane P	Positive control.
Lane N	Negative control.
LaneM	RBC 100 bp DNA Ladder Marker.



Figure 2. Nested PCR results of HCV at different concentrations. Viral nucleic acids were purified from samples containing different amount of HCV using MagCore® Viral Nucleic Acid Extraction Kit.

MagCore® Automated Nucleic Acid Kits Specification

MagCore®Super/HF16Plus/PlusII/EDA MagCore® HF16 / Compact / HF48 Cat No. Cat No. Cat No. Cat No. Cartridge Code Running Time Running Time 36 preps 96ргерз 36 preps 96ргерѕ

202

MagCore® Viral Nucleic Acid Extraction Kit (Low PCR Inhibition)

For 200 and 400 µl sample volumes.
Contents: Pre-Filled Cartridges, Proteinase K, PK Storage Buffer, Carrier RNA, Rnase Free Water, Disposable Tip & Holder Sets, Sample Tubes, Elution Tubes Shelf life: 18 months

62 min (sample volume: 200 µl) 57 min (sample volume:200 µl) MVN-400-04 73 min (sample volume:400 µl) 66 min(samplevolume:400 µl)

MagCore®Viral Nucleic Acid Large Volume Extraction Kit (1.2 ml)

For 1200 µl sample volumes.

Contents: Pre-Filled Cartridges, Proteinase K, PK Storage Buffer, Carrier RNA, Rnase Free Water, Disposable Tip & Holder Sets, Sample Tubes, Elution Tubes Shelf life: 18 months

80 min (sample volume:1200 µl) MVN-1200 MVN-1200 73 min (sample volume:1200 µl)

Enzyme Selection Guide

Product	Contents	Cat.No.
Proteinase K Set	11 mg Proteinase K,1.25 ml PK Storage Buffer	PK011
Carrier RNA Set	1 mg Carrier RNA, RNase Free Water	CR001











For Extraction of Viral DNA/RNA From Serum, Plasma, Swabs and Cell-Free Body Fluids by Using MagCore® System

MagCore® Viral Nucleic Acid Kits are designed to isolate viral DNA and/or RNA from plasma, serum, swabs and cell-free body fluids. All plastic consumables included in the kits are DNase/RNase-free to minimize the risk of cross contamination. Our patented magnetic beads will bind to the short DNA fragments with high affinity ensuring that all important genetic information is collected efficiently making your daily isolation routine easier and faster. This kit also features Internal Control selection in its protocol, being compatible with any viral detection kit.

203

MagCore® Viral Nucleic Acid Extraction Kit (High Sensitivity)

Magnetic beads
Separation Well 2

Magnetic beads
Separation Well 1

Empty
Lysis Buffer 1000µl

Binding Buffer 1000µl

Binding Buffer 1000µl

Wash 1 Buffer 1000µl

Wash 2 Buffer 1000µl

DEPCWater 1000µl

Wash 2 Buffer 1000µl

Wash 2 Buffer 1000µl

Wash 2 Buffer 1000µl

Wash 2 Buffer 1000µl

Features

- 1. Higher analytical sensitivity.
- 2. Consistent and reproducible results.
- 3. High performance of purified DNA/RNA in downstream applications such as aPCR.
- 4. Cartridges are pre-filled and sealed to prevent contamination.
- 5. Provides the option of Internal Control selection.


Applications

MagCore® Viral Nucleic Acid Extraction Kit allows automated processing of multiple sample types in the same run, including:

- 1. Cell-Free Body Fluids
- 2. Plasma and Serum
- 3. CSF
- 4. Urine
- 5. Swabs




Viral Nucleic Acid Extraction Performance Analysis

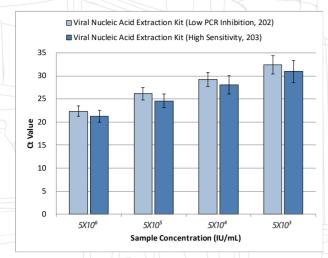

MagCore 2 203 shows better performance on real-time PCR with approximately $1\sim$ 2CtValueless ($n\geq$ 3)

Comparison of real-time PCR results of nucleic acid extraction by MagCore®, Brand Q and Brand R

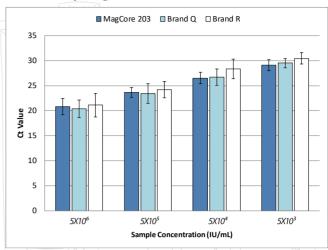

Purified viral nucleic acid by MagCore® is ideal for the downstream real-time PCR. Viral samples processed by MagCore®

Internal Control (IC) Selection

7_				1 111 111				
				Sample Concentration (IU/mL)—HBV Signal/IC Signal				
	No.	Kit	Signal	5x10 ⁶	5x10⁵	5x10 ⁴	5x10³	NTC
	1		HBV	22.06	25.39	28.58	30.53	
1		44 6 9202	+IC	29.99	31.54	31.17	31.55	4-
	2	MagCore®203	HBV	23.82	26.51	29.78	33.17	7-
			+IC	31.19	30.79	30.51	32.73	

MagCore®203 program provides internal control (IC) selection, and the real-time PCR analyzed by CE IVD HBV Quantification kit shows IC signal within criteria (Ct value 30±3). This confirms success in MagCore®203 viral extraction step and low co-purification of PCR inhibitors that may cause false amplification patterns.

MagCore® Automated Nucleic Acid Kits Specification

MagCore®Super/HF16Plus/PlusII/EDA						MagCore® HF16 / Compact / HF48			
Cautidas Cada	Cat No.		Cat No.	D. vereire e Tierre	Cat No.		CatNo.	Di versi in a Tinana	
Cartridge Code	36 preps		96preps	Running Time	36 preps		96preps	Running Time	

203 MagCore® Viral Nucleic Acid Extraction Kit (High Sensitivity)

For 200 and 400 µl sample volumes.

Contents: Pre-filled Cartridges, Proteinase K, PK Storage Buffer, Carrier RNA, RNase Free Water, Disposable Tip & Holder Sets, Sample Tubes, Elution Tubes.

Shelf life: 12 months

MVN400-05		MVN400-06	57 Mins/With IC Selection: 58 Mins (sample volume: 200 µl) 66 Mins/With IC Selection: 67 Mins (sample volume: 400 µl)	MVN400-05		MVN400-06	56 Mins/With IC Selection: 57 Mins (sample volume: 200 µl) 72 Mins/With IC Selection: 73 Mins (sample volume: 400 µl)
-----------	--	-----------	--	-----------	--	-----------	--

Enzyme Selection Guide

Product	Contents	Cat.No.
Proteinase K Set	11 mg Proteinase K, 1.25 ml PK Storage Buffer	PK011
Carrier RNA Set	1 mg Carrier RNA, RNase Free Water	CR001













For extraction of plant DNA from plant tissues

MagCore® Genomic DNA Plant Kit is designed to extract DNA from plant tissues and cells using MagCore® automated extraction systems. The provided Filter Column Set can filtrate hard tissue samples to prevent tissue residues from clogging the pipette syringe. The kit contains all required reagents and labware for automated extraction using magnetic-particle technology. Reagents are supplied and pre-filled in cartridges, which can be easily loaded into the MagCore® instrument.

301

MagCore® Genomic DNA Plant Kit



Features

- 1. Lysis buffer can degrade large amounts of polysaccharides present on plant cell walls
- 2. Cartridges are pre-filled and sealed to prevent contamination.
- 3. No phenol or chloroform extraction and alchol precipitation.
- 4. Efficient removal of contaminants and inhibitors.

Applications

The MagCore® Genomic DNA Plant Kit enables DNA purification from various samples types, including:

- 1. Plant cells
- 2. Plant tissues

High quality DNA available for downstream applications, including:

- 1. PCR and real-time PCR
- 2. Next Generation Sequencing (NGS)





Yields of nucleic acid purified from various sources by MagCore® Genomic DNA Plant Kit

Sample ID	Position	Yield (ng/μl)				
Citrus nobilisLour.	Leaf	36.5				
Carica papaya	Leaf	49.2				
Lycopersicon esculentum	Leaf	71.9				
Ctrullus lanatus	Leaf	23.4				
Anisogonium esculentum	Leaf	240.3				
Trachycarpus fortunei	Leaf	154.7				
Mentha piperita	Leaf	54.6				
Anthoceros punctatus	Leaf	38.9				
Calix babylonica	Leaf	80.7				
Prnus campanulata	Leaf	35.7				
Lxeris chinensis	Leaf	161.4				
Capsicum annuum	Leaf	13.9				
Vigna radiata	Seed	7.0				
Medicago sativa	Leaf	21.0				
Vigna angularis	Seed	5.8				
Agaricus bisporus	Whole	14.3				

MagCore® Automated Nucleic Acid Kits Specification

	N	lagCore®Super/HF1	MagCore® HF16 / Compact / HF48				
Cartridge Code	CatNo.	CatNo.	Punning Timo	CatNo.	CatNo.	Dunning Time	
	36 preps	96preps	RunningTime	36 preps	96preps	Running Time	

MagCore® Genomic DNA Plant Kit
For 50-100 mg fresh tissues or 5-20mg dried plant tissues
Contents: Pre-Filled Cartridges, Rnase A, GP1 Buffer, GP2 Buffer, Filter Column Sets, Disposable Tip & Holder Sets, Sample Tubes, Elution Tubes
Shelf life: 18 months

MCD 01	MGP-01	MGP-02	Time:33 min (sample volume:400 µl)	MGP-01		MCD 02	33min(sample volume:400 µl)
IVIGP-UI			*optical detection is not provided	MGP-01		MGP-02	(Applicable models : HF16, Compact)

Enzyme Selection Guide

Product	Contents	Cat.No.
Proteinase K Set	11 mg Proteinase K, 1.25 ml PK Storage Buffer	PK011
	50 µl RNase A (50mg/ml)	RN050
RNase A	130μlRNaseA (50mg/ml)	RN130









RBC Bioscience Corp. www.rbcbioscience.com info@rbcbioscience.com





For extraction of genomic DNA from tissues, swabs, FFPE, stool, feed/soil and forensic samples

40 Genomic DNA Tissue Kit is designed for purification of total DNA (including genomic, mitochondrial and viral DNA) from a variety of animal tissues or cells by using MagCore® auto-extraction instrument. The provided Filter Column can filtrate hard tissue sample or swab sample to prevent tissue residues to obstruct pipette tip during the process of MagCore®. The method uses pre-filled cartridges based on cellulose coated magnetic bead technology.

401

401 MagCore® Genomic DNA Tissue Kit



Features

- 1. Suitable for a wide variety of sample types, including forensic samples
- 2. Consistent and reproducible results
- 3. Purified DNA of approximately 20-30 kb in length is suitable for PCR or other enzymatic reactions.

Applications

MagCore®Genomic DNA Tissue kit allows processing multiple sample types in the same run, including:

- 1. Ānimal Tissues: Solid animal tissue, Stool samples
- 2 Swah
- 3. Feed and soil samples
- 4. Formalin-fixed, Paraffin-embedded (FFPE) tissues samples (manual pretreatment)
- 5. Cultured Yeast samples
- 6. Forensics samples



DNA yield gained from 10 mg of mice tissue samples and 0.25 cm tail

Mice tissue	Yield (ng/ul)
Liver	17.5
Spleen	17.5
Brain	10
Lung	12.5
Kidney	32.5
Tail	7.5

Table 1. Genomic DNA was purified from 10mg of mice tissue samples and 0.25cm tail.

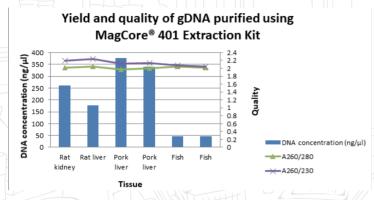


Figure 1. Genomic DNA from 10mg of various animal tissues were extracted by the MagCore®401 Genomic DNA Tissue Kit. DNA was eluted in 60 µL and quantitated using Nanodrop n-1000.

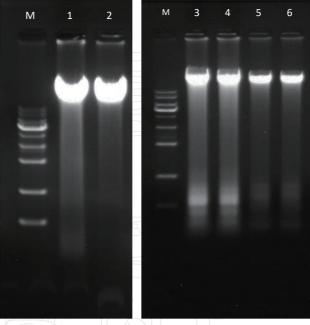


Figure 2. Agarose gel electrophoresis of the various animal tissue DNA was isolated from MagCore® automated extraction.; Lane M: RBC 1 kb DNA ladder Marker; Lane 1:Ratkidney;Lane 2:Ratliver;Lane 3-4:Porkliver;Lane 5-6:Fish tissue.

MagCore® Automated Nucleic Acid Kits Specification

MagCore®Super/HF16Plus/PlusII/EDA					MagCore® HF16 / Compact / HF48		
Cartridge Code	CatNo.	Cat No.	RunningTime	Cat No.		CatNo.	RunningTime
	36 preps	96preps		36 preps		96preps	

401

MagCore®Genomic DNA Tissue Kit

For tissue and forensic samples (up to 40mg)
Contents: Pre-Filled Cartridges, Filter Column Sets, GT Buffer, Proteinase K, PK Storage Buffer, Disposable Tip & Holder Sets, Sample Tubes, Elution Tubes Shelf life: 18 months

MGT-01	MGT-02	33 min (samplevolume:400 µl) * optical detection is not provided	MGT-01	MGT-02	33 min (sample volume:400 μl)

Enzyme Selection Guide

Product	Contents	Cat.No.
Proteinase K Set	11 mg Proteinase K, 1.25 ml PK Storage Buffer	PK011
DM 4	50 µl RNase A (50mg/ml)	RN050
RNase A	130 µIRNase A (50mg/ml)	RN130









RBC Bioscience Corp. www.rbcbioscience.com info@rbcbioscience.com





For Automated Extraction of Genomic DNA from Formalin-fixed Paraffin-embedded (FFPE) Tissue

MagCore® Genomic DNA FFPE One-Step Kit is designed to purify total DNA (including genomic, mitochondrial, and viral DNA) from Formalin-fixed Paraffin-embedded tissues via MagCore® Automated Extraction instruments. Our program features One-Step Heating, which automatically melts paraffin and lyses cells at the same time. No harmful reagent such as xylene is involved in the deparaffination process. Two protocols are outlined for different tissues sizes: 2-hour setup for smaller tissues and 16-hour setup for larger tissues.

405

MagCore® Genomic DNA FFPE One-Step Kit

Magnetic beads
Separation Well 2

Magnetic beads
Separation Well 1

GT buffer 500µl
Lysis Buffer 1000µl
Binding Buffer 1000µl
Beads Mixture 500µl
Wash 1 Buffer 1000µl
Wash 2 Buffer 1000µl
Wash 2 Buffer 1000µl
Elution Buffer 1000µl
Elution Buffer 1000µl
Empty
Empty
Empty
Heat block Well 1

Features

- 1. Can load FFPE tissue samples with a surface area of up to 300 mm2 and up to 5 µm thick.
- A full package for DNA to be purified from FFPE samples. Complete automatic procedure from deparaffination to elution.
- 3. Stable deparaffination by Sula oil and reproducible results; no xylene involved.
- 4. Thermostavle cap is specially designed to cover the reaction well for an optimal condition where the lysis buffer and the FFPE sample can mix properly during the processing period.

Applications

- 1. PCR and real-time quantitative PCR
- 2. Southern Blot
- 3. Next Generation Sequencing
- 4. Genotyping



2-hr heating protocol works out


There are 2-hour heating and 16-hour heating protocols. It is suggested for a better performance that the 2-hour heating program is for small tissues and the 16-hour program is for large tissues.


Genomic DNA with high performance

The DNA extracted by MagCore® Genomic DNA FFPE One-Step Kit is high yield and suitable for downstream PCR and real-time PCR.

A. DNA concentration


Whether samples were treated with either a 2-hour or 16-hour heating, yields of DNA extracted by MagCore® are mostly higher than the yields of DNA extracted by the competitor.

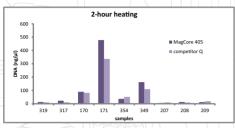

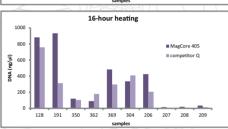

B. The downstream PCR

PCR products of DNA purified by MagCore® and the competitor brand are confirmed by 1.5% agarose gelelectrophoresis.

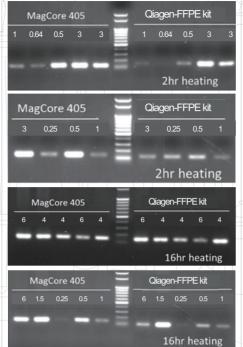



DNA yield gained from 10 mg of mice tissue samples and 0.25 cm tail

OLED CITTOM	_//		
Sample ID			Yield (ng/ul)
Liver			17.5
Spleen			17.5
Brain			10
Lung			12.5
Kidney			32.5
Tail			7.5

C. Purified DNA ideal for the downstream real-time PCR

Whether samples experience a 2-hour or 16-hour heating, the CT values of samples processed by MagCore® are mostly lower than the CT values of samples processed by the competitor.

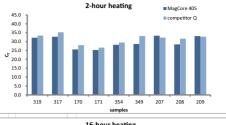

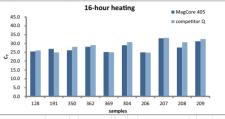

D. Tissue sample sizes

It is shown below that different sizes of tissue samples were processed by a 2-hour or 16-hour heating program.

2-hour heating								
	Sample	Size(cm²)						
1	319	1*1						
2	317	0.8*0.8						
3	170	0.5*1						
4	171	1.5*1						
5	354	1.5*2						
6	349	1.5*2						
7	207	0.5*0.5						
8	208	0.5*1						
9	209	0.5*1						

	16-hourheating						
	Sample	Size(cm²)					
1	128	2*3					
2	191	<i>2</i> *2					
3	350	2*2					
4	362	2*3					
5	369	<i>2</i> *2					
6	304	2*3					
7	206	1.5*1					
8	207	0.5*0.5					
9	208	0.5*1					
10	209	1*1					

MagCore® Automated Nucleic Acid Kits Specification

_	· · · · · · · · · · · · · · · · · · ·							
		MagCore®Super/HF10	MagCore® HF16 / Compact / HF48					
Cartridge Code	Cat No.	CatNo.	RunningTime	Cat No.	CatNo.	Dunaina Tima		
	36 preps	96preps		36 preps	96preps	RunningTime		

405

MagCore® Genomic DNA FFPE One-Step Kit

For FFPE tissue samples.

Contents: Thermostable Cap, Sula Oil, Pre-Filled Cartridges, Filter Column Sets, GT Buffer, Proteinase K, PK Storage Buffer, Disposable Tip & Holder Sets, Sample Tubes, Elution Tubes Shelf life: 18 months

MGF-01		MGF-03	33min(samplevolume:400 µl)	MGF-01		MGF-03	175 min(2-hour heating)-Standard 1012 min(16-hour heating)-High Yield
--------	--	--------	----------------------------	--------	--	--------	--

Enzyme Selection Guide

	Product	Cat.No.	
	Proteinase K Set	11 mg Proteinase K,1.25 ml PK Storage Buffer	PK011
211	50µl RNaseA (50mg/ml)	RN050	
	130 µI RNase A (50mg/ml)	RN130	









RBC Bioscience Corp. www.rbcbioscience.com info@rbcbioscience.com





For extraction of genomic DNA from forensic samples

MagCore® Genomic DNA Forensic Direct Kit is designed for purifying total DNA from forensic samples, such as dried blood spots, cigarette butts, cartilage, hair roots, seminal stains, and chorionic villus, using MagCore® auto-extraction instruments. Its unique feature is RBC patented technology that allows to isolate DNA automatically from solid samples without any pretreatment.


406

MagCore® Forensic DNA Direct Kit




Applications

It uses magnetic-particle technology to purify genomic DNA from forensic samples. The purified genomic DNA can be directly used for downstream applications such as STR, PCR and real-time PCR.

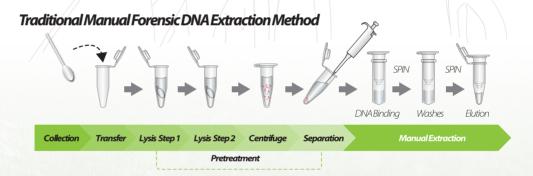


Automated Lysis and Separation

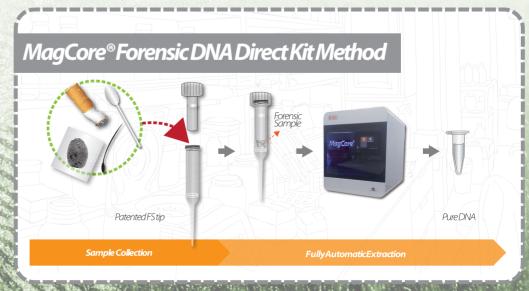



Unique Pipette Tip Design

- Airtight assurance (hermetic seal)
- Automatic separation of the solid debris

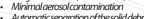

Compare the results between 406 and Q kit by qPCR



Other casework samples



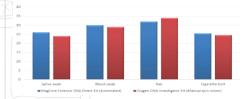


Figure: Extraction of simulated casework samples. This test was analyzed by real-time PCR with GAPDH gene primer. Salivs swab: 10 ml of saliva was applied (n=4), Blood swab: 1 ml of blood was applied and allowed to dry (n=4), Hair: 1 hair each from donor (n=4), Cigarette butt: A quarter of a filter paper (n=4).



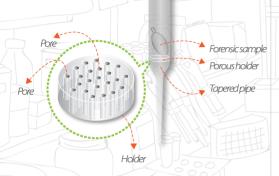

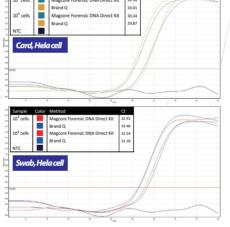

MagCore® Automated Nucleic Acid Kits Specification

Tip collars

Top filter

Designed hold

MagCore®Super/HF16Plus/PlusII/EDA

Cat No. Cartridge Code Running Time 72preps

406

MagCore®Forensic DNA Direct Kit

For extracting genomic DNA from forensic samples
Contents: Pre-Filled Cartridges (Including Proteinase K), Disposable Tip & Holder Sets, Elution Tubes, FS Tip, 200 ml SP Tip

MFC-03 120min

Enzyme Selection Guide

Shelf life: 12months

Product	Contents	Cat.No.
Proteinase K Set	11 mg Proteinase K, 1.25 ml PK Storage Buffer	PK011
201/	50 µl RNase A (50mg/ml)	RN050
	130µIRNaseA (50mg/ml)	RN130









RBC Bioscience Corp. www.rbcbioscience.com info@rbcbioscience.com





For extraction of genomic DNA from bacteria

MagCore®Genomic DNA Bacterial Kit is designed to extract genomic DNA from Gram positive and Gram negative bacteria. The Kit contains all required reagents and labware for automated extraction using magnetic-particle technology. Reagents are supplied and prefilled in cartridges, which can be easily loaded into the MagCore® instrument.

502

MagCore® Genomic DNA Bacterial kit

Magnetic beads
Separation Well 2

Magnetic beads
Separation Well 1

Empty
Lysis Buffer 1000µl
Binding Buffer 1000µl
Beads Mixture 500µl
Wash 1 Buffer 1000µl
Beads Mixture 1000µl
Empty
Empty
Empty
Empty
Empty
Empty


Features

- 1. High performance of purified DNA in downstream applications such as gPCR.
- 2. Cartridges are pre-filled and sealed to prevent contamination.
- 3. No phenolor chloroform involved.

Applications


MagCore® Genomic DNA Bacterial Kit is designed for automated DNA extraction from various sample types, including:

- 1. Cell cultures in suspension (up to 5x10° cells)
- 2. Biological fluids
- 3. Cell cultures on plate
- 4. Gram-positive and Gram-negative bacteria

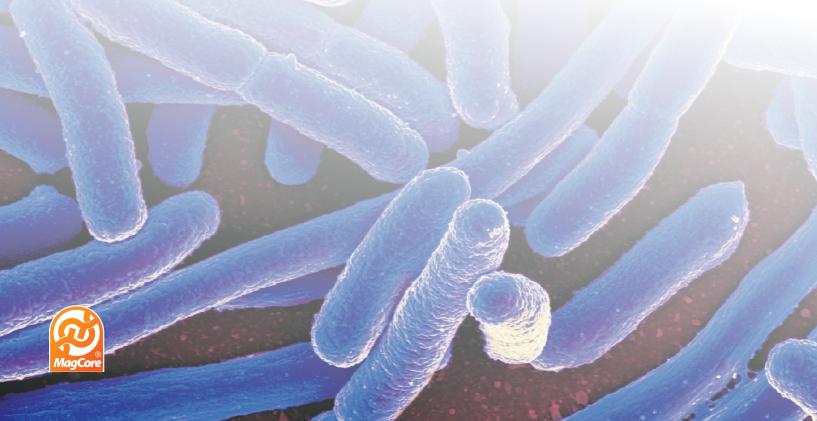


High performance of purified bacterial DNA in real-time PCR

Well	Туре	Target	Ct	Interpretation	Label
D2	Unknow	H. pylori	30.70	Positive	Sample A
D3	Unknow	H. pylori	30.52	Positive	Sample A
D4	Unknow	H. pylori	No Ct	Negative	Negative Ctrl
D5	Unknow	H. pylori	30.59	Positive	Sample A
E2	Unknow	H. pylori	26.73	Positive	Sample B
E3	Unknow	H. pylori	26.28	Positive	Sample B
E4	Unknow	H. pylori	26.40	Positive	Sample B

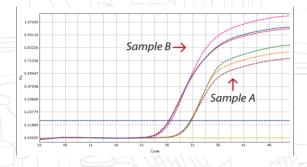

Consistent Yield, Purity, and Reproducibility in MagCore® Bacteria Genomic DNA Extraction



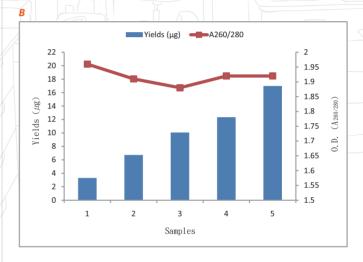

MagCore® Automated Nucleic Acid Kits Specification

MagCore®Super/HF16Plus/PlusII/EDA MagCore® HF16 / Compact / HF48 Cat No. Cat No. Cat No. CatNo. Cartridge Code Running Time Running Time 36 preps 96preps 96preps 36 preps

502

MagCore®Genomic DNA Bacterial kit

For 1200µl sample volumes. Contents: Pre-Filled Cartridges, Lysozyme Reaction Buffer, Proteinase K, PK Storage Buffer, Rnase A, Disposable Tip & Holder Sets, Sample Tubes, Elution Tubes Shelf life: 18 months

MBB-01		MBB-02	39 min (sample volume:200 µl) * optical detection is not provided	MBB-01		MBB-02	44min(samplevolume:200μl)
--------	--	--------	--	--------	--	--------	---------------------------

Enzyme Selection Guide

Product	Cat.No.	
Proteinase K Set	PK011	
nnaseA	50µl RNaseA (50mg/ml)	RN050
	130 µl RNase A (50mg/ml)	RN130









RBC Bioscience Corp. www.rbcbioscience.com info@rbcbioscience.com





For extraction of microbial and host genomic DNA from stool samples

MagCore® Gut Microbiome DNA Kit is designed specifically for isolating high-quality microbial and host genomic DNA from stool and gut samples. With a short pretreatment, DNA from tough-to-lyse Gram-positive and Gram-negative bacteria which inhabit fecal samples can be successfully isolated with efficient lysis and homogenized with Beads-Beating Tube. In addition, inhibitor substances (e.g. polysaccharides, protein, etc.) commonly found in stools are removed with one-step IR Buffer. The crude DNA is then precipitated on ice and resuspended to adjust the condition that is suitable for binding genomic DNA.

Extracted DNA is ready for downstream molecular-based applications including qPCR, 16S rRNA gene sequencing, etc.

504

MagCore® Gut Microbiome DNA Kit

Magnetic beads
Separation Well 1

Magnetic beads
Separation Well 1

Proteinose K50,J

Empty
Binding Buffer GM
15000,J

Wash 1Buffer 1000,J

Wash 2Buffer 1000,J

Bution Buffer 1000,J

Empty
Empty
Empty
Empty
Empty

Features

- Streamlined protocols for the isolation of inhibitor-free, PCR-quality DNA (up to 50 μg/prep) from microbes including Gram-positive and Gram-negative bacteria.
- 2. Hands-on time can be as little as 20 minutes.
- 3. Depletion of PCR inhibitors by precipitation using a novel one-step IR Buffer.
- 4. High-purity DNA is ready for common downstream applications such as qPCR and 16S rRNA gene sequencing.


Applications


The purified DNA can be directly used for downstream application such as:

- 1. qPCR
- 2. 16S rRNA gene Sequencing
- 3. Next-Generation Sequencing Applications




Stool DNA Extraction Efficiency

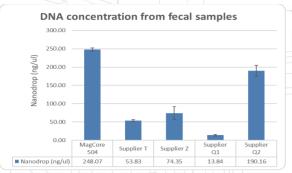

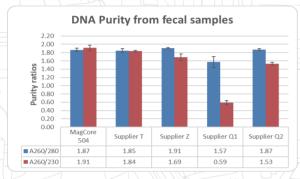


DNA was isolated from 0.2 g stool samples (in triplicates) with the MagCore Gut Microbiome DNA kit and supplier kits. (supplier T, supplier Z, supplier Q1, supplier Q2). DNA concentration and purity (A260/280, A260/230) are measured by Nanodrop spectrophotometer.

Gel electrophoresis

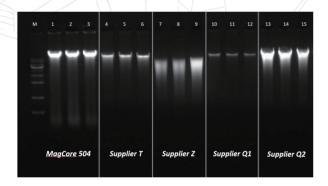


Analysis of Genomic DNA extracted from human stool samples using 1% gel electrophoresis. Performance of MagCore Gut Microbiome kit analyzed on agarose gel electrophoresis; Lane M: RBC 1 kb DNA ladder Marker.

16SrRNA gene qPCR

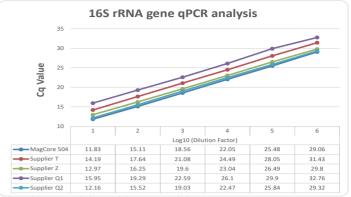


DNA extracted from stool samples using MagCore 504 kit and other competitor kits were evaluated by 10-fold serial dilution of 16S rRNA gene qPCR amplification. Non template control (NTC) was included as a reference. (amplicon size: 466bp).

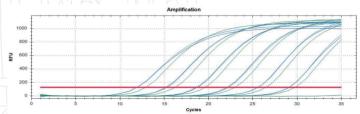

qPCR amplification curve of MagCore® 504, supplier Q2, supplier Z: MagCore 504 shows a low inhibition, similar amplification curve with supplier Q2, and superior Cq values to supplier Z.

Human fecal microbiota qPCR analysis

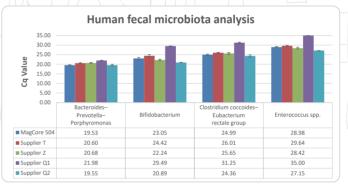

Phylum	lum Target organism				
Bacteroidetes	Bacteroides-Prevotella-Porphyromonas	G-			
Actinobacteria	Bifidobacterium	G+			
Actinobacteria	Clostridium coccoides-Eubacterium rectale group	G+			
Firmicutes	Enterococcus spp.	G+			

Human fecal microbiota qPCR analysis of four bacterial targets: Bacteroides-Prevotella-Porphyromonas, Bifidobacterium, Clostridium coccoides-Eubacterium rectale group, Enterococcus spp.. DNA from triplicate stool samples were isolated wih MagCore gut microbiome DNA kit and other suppliers.

MagCore® Automated Nucleic Acid Kits Specification

MagCore®Super/HF16Plus/PlusII/EDA						MagCore® HF16 / Compact / HF48			
Cartridge Code	Cat No.	Cat No.		RunningTime	Cat No.	CatNo.		Dunaning Time	
	24preps	72preps			24preps	72preps		RunningTime	

504

MagCore®GutMicrobiomeDNAKit

For 200±20 mg stool sample input.

Contents: Pre-filled cartridges, Beads-Beating Tube, GM-L1 Buffer, GM-L2 Buffer, GM-IR Buffer, GM-P1 Buffer, GM-P2 Buffer, GM-LY Buffer, Disposable tip & holder sets, Sample tubes, Elution tubes
Shelf life: 12 months

MGM-01	MGM-02	Mage	ore®Super–52min ore®Plus II-55min	MGM-01	MGM-02		
--------	--------	------	--------------------------------------	--------	--------	--	--











For extraction of total RNA from formalin-fixed paraffin-embedded (FFPE) tissue by using MagCore® System.

MagCore® Total RNA FFPE One-Step Kit is specially designed for total RNA purification from FFPE tissues by MagCore® instruments. It features the method, one-step heating, to melt paraffin without harmful reagents involved such as xylene or other organic solvents, and lyse tissues at the same time. The MagCore® Total RNA FFPE One-Step Kit System optimizes the lysis conditions to reverse the formalin fixation without the need for overnight digestion and retain both large and small RNAs. The program provides optional DNase I treatment to remove contaminated DNA.



MagCore® Total RNA FFPE One-Step Kit



Features


- 1. A full package for total RNA isolation from FFPE tissue samples. It starts from melting paraffin to the final RNA
- 2. Highly user-friendly protocol with minimal pretreatment required.
- 3. Sula óil (Deparaffinization solution): no xyléne or other organic solvents involved.
- 4. Thermostable cap is especially designed to cover the reaction well for an optimal condition where lysis buffer and the FFPE samples can mix properly during the long processing period.
- 5. The total processing time is within approximately 160 mins.

Applications

The purified RNA can be directly used for downstream application such as:

- 1. cDNA synthesis
- 2. real-time PCR
- 3 RT-PCR
- 4. Microarray
- 5. Next-Generation Sequencing

FFPERNA extraction efficiency

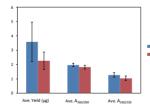


Rat tissue	Average Yields (µg)	Average A ₂₆₀ / ₂₈₀	Average A ₂₆₀ / ₂₃₀		
Liver	3.57±1.38	1.96 ±0.11	1.26±0.16		
Kidney	2.25±0.62	1.82±0.12	1.03 ±0.15		


Total RNA was purified from various FFPE rat tissues stored at -80°C for 3 months. RNA yield and quality from one (liver, kidney) 50 µm tissue section per sample was determined by Nanodrop n-1000.


1.89

2.17

MagCore®605 competitors

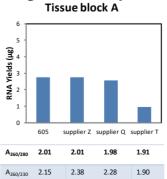

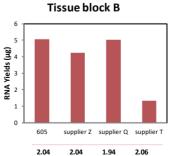

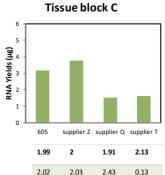


2.43


MagCore® 605 retains high RNA extraction efficiency in various rat liver tissue blocks compare to several FFPE RNA purification kit suppliers. The RNA yield and quality was quantified by Nanodrop'n -1000.


2.3

2.08

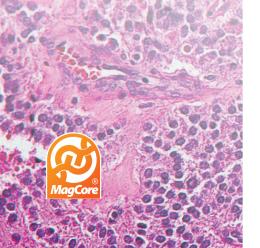

RNA integrity

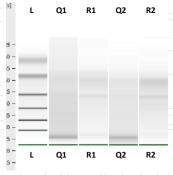

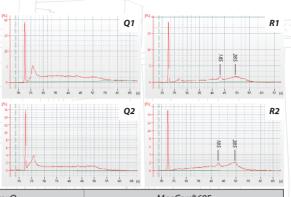


1 1-11							
		Competitor Q		MagCore® 605			
Tissue Block	Sample	RNA yields (ng/µl)	RIN	Sample	RNA yields (ng/μl)	RIN	
#1	Q1	97.37	2.2	R1	<i>77.</i> 9	3.2	
#2	Q2	<i>58.7</i>	2.2	R2	62.6	3.3	

RNA integrity assessment

Total RNA isolation from rat liver 50µm tissue sections of FFPE tissue blocks were measured on Agilent 2100 bioanalyzer. Isolated RNA from MagCore® 605 kit maintains its integrity (18S, 28S peaks).

MagCore® 605 competitors

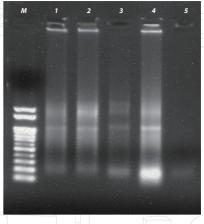


Brand	Sample	nanodrop (ng/µl)	A ₂₆₀ / ₂₈₀	A ₂₆₀ / ₂₃₀	Ct Value (GAPDH)
MagCore®	1	60.12	1.93	1.48	24.01
605	2	74.79	1.89	1.52	22.02
SupplierZ	3	31.21	1.96	1.49	26.61
Supplier Q	4	117.54	1.93	2.17	23.75
SupplierT	5	5.89	1.46	0.15	27.39

Comparison of total RNA extraction efficiency by using various FFPE RNA extraction kit suppliers. Total RNA was purified from one rat liver 50 µm tissue section of a single FFPE tissue block for each sample. The RNA yield and quality was quantified by Nanodropn-1000.

RT-PCR performance



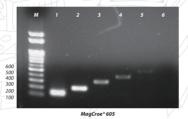





Evaluation of RNA performance in RT-PCR

Five amplicons of different length (118, 206, 312, 400, 503, 613 bp amplicon size) of the gene Rpl4 performed using cDNA derived from FFPE rat liver samples purified by various suppliers.

MagCore®Automated Nucleic Acid Kits Specification

MagCore®Super/HF16Plus/PlusII/EDA						Мо	agCore® HF16	:/Compact	
CartridasCodo	Cat No.	CatNo.		Running Time	Cat No.	CatNo.		Di vanina Timo	
Cartridge Code	24preps	72preps			24preps	72preps		RunningTime	



MagCore®TotalRNAFFPEOne-stepKit

For extraction of total RNA from formalin-fixed paraffin-embedded (FFPE) tissue by using MagCore® System. Contents: Pre-Filled Cartridges, Disposable Tip & Holder Sets, Elution Tubes Shelf life: 12 months

MRF-01 MRF-03	140min (without DNase I treatment) 158min (with DNase I treatment)	MRF-01	MRF-03		147min (without DNase I treatment) 165min(with DNase I treatment)
---------------	---	--------	--------	--	--

Enzyme Selection Guide

Product	Contents	Cat.No.
BW 45	For 36 reactions RNase-Free DNase I (Lyophilized):1500 Kunitz units x 1 vial, 1 ml RNase-Free Water x 1, 15 ml DNase I Reaction Buffe	DN036
DNase I Set	For 96 reactions RNase-Free DNase I (Lyophilized): 1500 Kunitz units x 2 vial, 1 ml RNase-Free Water x 2, 30 ml DNase I Reaction Buffer	DN096














For purification of miRNA from serum and plasma


MagCore® Plasma miRNA Extraction kit is designed for purification of high purity microRNA from 400µl serum and plasma using MagCore® automated extraction instrument. All the kit components of plastic consumables are DNase/RNase-free pretreated, to eliminate the possibility of cross contamination.




High recovery of microRNA

 $Extraction \ results of microRNA isolation from 400 \mu l plasma from 3 \ different donors \ with an elution volume of 60 \mu l using MagCore Plasma \ miRNA Extraction \ kit (620) \ or \ Brand \ Q \ kit.$

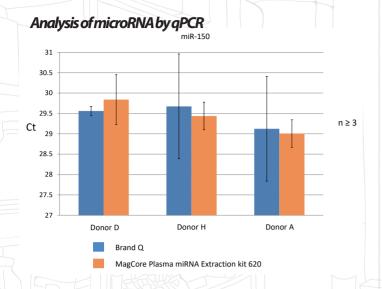

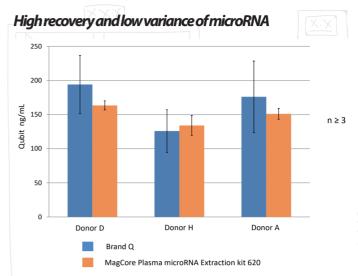

Detection of microRNAs

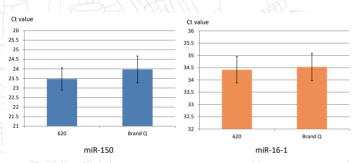

micrioRNA was isolated from 400 µl plasma MagCore Plasma miRNA Extraction kit (620) or Brand Qkit. miR-150 and miR-16-1 was detected by Real-Time PCR.



MagCore®Automated Nucleic Acid Kits Specification

				,					
MagCore® Super/HF16 Plus/Plus II/EDA						MagCore® HF16/Compact			
CartridasCodo	CatNo.	CatNo.		Di vanina Timo	CatNo.	CatNo.	RunningTime		
CartridgeCode	24 preps	72preps		Running Time	24 preps	72preps	runing inne		
620	For 400 µl sa	mple volumes -Filled Cartridge		ktraction Kit Buffer(10 ml), Disposable Pipet Tip plus	Holder Sets, Samp	ple Tubes, Elution Tub	ne		
	MMR-01	MMR-03		40 min (without DNase I treatment)	MMR-01	MMR-03	34 min (without DNase I treatment)		

60 min (with DNase I treatment)









57 min (with DNase I treatment)





For total RNA extraction from cultured cells, whole blood and tissues

MagCore® triXact RNA Kit is specially designed for total RNA purification from the three most common sample types in diagnostics, research, and forensics: up to 1x10⁶ cultured cells, a variety of tissues, or whole blood samples. The program provides optional DNase I treatment to remove residual DNA and extract high quality DNA-free RNA. Make use of its oustanding extraction performance, effiency, plus easy-to-follow protocols to simplify your every day extraction!



MagCore®triXact RNA Kit

Separation Well 2
Separation Well 2
Separation Well 1
Empty
Lysis Buffer 1000µl
Binding Buffer 1000µl
Beads Mixture 1000µl
Wash 1 Buffer 1000µl
Wash 2 Buffer 1000µl
DEPCH,O 1000µl
DEPCH,O 1000µl
DEPCH,O 1000µl
DEPCH,O 1000µl
Heart block Well 2
Heart block Well 2

Features

- 1. Excellent RNA Yield and Purity shown in both A_{260}/A_{280} ratio and RIN value
- 2. Highly efficienct and user-friendly protocol with minimal pretreatment required
- 3. Capable of isolating RNA from 3 types of samples: Cells, Whole blood, Tissues
- 4. Optional DNase I treatment to remove residual DNA

Applications

Purified RNA ready to use in downstream applications such as:

- 1. PCR
- 2. Real-time PCR
- 3. Microarray target preparation
- 4. Northern blotting
- 5. NGS analysis



Total RNA from 1 x 10⁶ Hela Cell Extract

No.	Kit	nanodrop ng/μl	260/280	260/230
1		207.15	2.08	1.86
2	Brand Q	182.85	2.07	1.78
3		174.90	2.11	1.63
4	Brand Z	171.52	2.03	2.24
5		210.45	2.02	2.22
6		164.76	2.05	2.17
7	MagCore*	264.69	2.04	2.11
8	triXact RNA Kit	250.61	2.02	2.17
9		237.44	2.03	2.17
10	Brand M	224.35	2.11	2.26
11		252.69	2.10	2.30
12		222.45	2.08	2.18



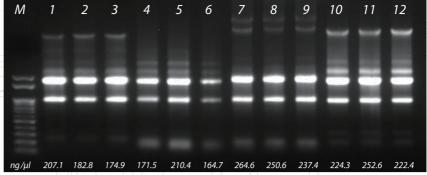


Fig 1. Gel Electrophoresis of RNA from 1 x 106 Hela Cell Extract

ng/μl		Concentr	ation	nanodrop
20.00				
00.00			I	
80.00				
60.00				
40.00				
20.00	I	I		
0.00 +	Brand Q	Brand Z	MagCore® triXact RNA	Brand M

Fig 3. Concentration comparison of RNA from 400µl Human Whole Blood

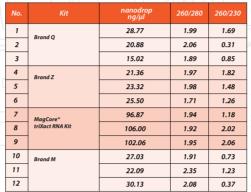


Table 2. Analysis of Whole Blood RNA isolated by MagCore $^{\!\circ}$ triXact RNA Kit and others

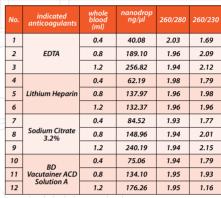


Table 3. Extraction efficiency of RNA from Human Whole Blood with various anticoagulants

Total RNA from 30 mg Fish Tissue

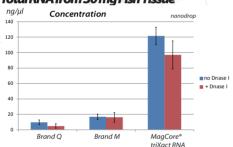


Fig 4. Concentration comparison of RNA from 30 mg Fish Tissue with/without DN as el treatment

Specification

MagCore®Automated Nucleic Acid Kits Specification

MagCore®Super/HF16Plus/PlusII/EDA

MaaCore®	HF16 /	Compact
MagCore®	HF 16 /	Compaci

Running Time

Cartridge Code

24preps	72preps	Running Time	24preps	72preps
Cat No.	Cat No.		Cat No.	CatNo.



For total RNA extraction from cultured cells, human whole blood and animal tissues

ror total in we know that the cens, in a man whole blood and a minial issues Contents: Pre-Filled Cartridges, RB Buffer, RBC Lysis Buffer, Filter column Set, Disposable Tip & Holder Sets, Sample Tubes, Elution Tubes Shelf life: 12 months

MRX-01	MRX-03		48min (without DNasel treatment; starting volume:400 µl) 81 min (with DNasel treatment; starting volume:400 µl) * optical detection is not recommended	MRX-01	MRX-03		S8min (without DNasel treatment; starting volume.400 µl) 73min (with DNasel treatment; starting volume.400 µl) (Applicable models: HF16, Compact.)
--------	--------	--	--	--------	--------	--	--

Enzyme Selection Guide

Product	Contents	Cat.No.
Proteinase K Set	11 mg Proteinase K, 1.25 ml PK Storage Buffer	PK011
	For 36 reactions RNase-Free DNase I (Lyophilized):1500 Kunitz units x 1 vial, 1 ml RNase-Free Water x 1, 15 ml DNase I Reaction Buffe	DN036
DNase1Si	For 96 reactions RNase-Free DNase I (Lyophillized): 1500 Kunitz units x 2 vial 1 ml RNase-Free Water x2, 30 ml DNase I Reaction Buffer	DN096













For extraction of poly(A)-mRNA from human whole blood, culture animal cell, total RNA

MagCore® mRNA One-Step Kit is specially designed for poly(A)-mRNA purification from human whole blood/culture animal cell/total RNA. The program offers optional DNase I treatment to remove residual DNA from contaminating the results. High quality DNA-free mRNA can be extracted using this kit along with our RNase-free DNase I. The isolation protocol relies on base pairing between the polyA residues at the 3' end of most eukaryote mRNA, and the oligo (dT)25 residues covalently coupled to the surface of the beads. Other RNA species lacking a polyA tail will not hybridize with the beads and are readily washed away.



MagCore®mRNA One-Step Kit



Features

- 1. Exclude the RNA species which lacks a polyA tail
- 2. Extract High-quality DNA-free mRNA by combining with our RNase-free DNase I
- 3. Optional DNase I treatment to remove residual DNA

Applications

Purified poly(A)-mRNA can be directly used for downstream application such as:


- 1. Real-time RT-PCR
- 2. RT-PCR
- 3. cDNA synthesis






Blood mRNA extraction efficiency

			400ul pre-treatment Blood					Elution
	Kit	without DNase		with DNase				
		Magcore 639			Magcore 639			buffer
mRNA 100nt	Average	22.10	21.53	22.12	22.32	21.83	25.05	N/A
	SD	0.14	0.26	0.26	0.21	0.48	0.08	
	CV	0.6%	1.2%	1.2%	0.9%	2.2%	0.3%	
mRNA 500nt	Average	22.86	22.05	22.95	23.10	22.54	27.06	N/A
	SD	0.33	0.89	0.43	0.24	0.53	1.58	
	CV	1.3%	4.0%	1.9%	1.0%	2.3%	5.8%	
DNA 100bp	Average	30.10	36.41	27.47	38.72	36.86	38.21	N/A
	SD	0.91	0.36	0.48	1.43		0.23	
	CV	3.0%	1.0%	1.7%	3.7%	-	0.6%	

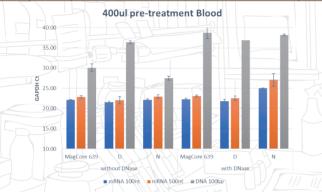


Cultured cell mRNA extraction efficiency

		1*10^6THP-1 cell						
	Kit	with	out DNase		with DNase			Elution
		Magcore 639			Magcore 639			buffer
	Average	17.18	17.37	16.48	17.1	17.18	17.17	N/A
mRNA 100nt	SD	0.62	0.35	0.02	0.2	0.18	0.19	
TOOTE	cv	3.60%	2.00%	0.10%	1.20%	1.00%	1.10%	
	Average	19.59	19.46	19.46	20	19.48	19.4	N/A
mRNA 500nt	SD	0.34	0.54	0.29	0.08	0.18	0.21	
Soork	cv	1.70%	2.70%	1.50%	0.40%	0.90%	1.10%	
	Average	32.15	34.85	26.48	36.96	37.8	40	
DNA 100bp	SD	0.5	1.56	0.42				N/A
	cv	1.50%	4.50%	1.60%	-	-	-	

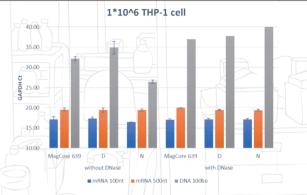

Efficient removal of ribosomal RNA

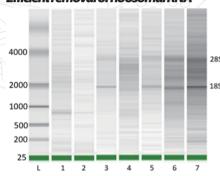

- L: Ladder
- 1: MagCore 639 mRNA from whole blood
- 2: MagCore 639 mRNA from pre-treatment blood
- 3: Dbrand kit mRNA from pre-treatment blood 4: Nbrand kit mRNA from pre-treatment blood
- 5: MagCore 639 mRNA from THP-1 cell
- **6:** Dbrand kit mRNA from THP-1 cell
- 7: NbrandkitmRNA from THP-1 cell

	Kitbrand	Sample	Automatic/Manual	rRNA contamination (%)	
1	Magcore639	Wholeblood	Automatic	0	
2	Magcore 639		Automatic	0	
3	D	Pre-treatment blood	Manual	6	
4	N		ivianuai	0	
5	Magcore 639		Automatic	2.6	
6	D	THP-1 cell	Manual	6.9	
7	N		ivianuai	2.8	

MagCore® Automated Nucleic Acid Kits Specification

MagCore®Super/Plus II/Plus II Dx/EDA

Cartridge Code

				1	 	
Cat No.	Cat No.		Running Time			
24preps	72preps					



MagCore®mRNA One-Step Kit

For extraction of poly(A)-mRNA from human whole blood, culture animal cell, total RNA
Contents: Pre-Filled Cartridges, mRNA beads Mixture (13mL), Pipet Tip plus Holder Set, Sample Tube, Elution Tube, Lysis buffer (mRNA), RBC lysis buffer (200 ml)
Shelf life: 12 months

MRM-01 MRM-03 S2min (without DNase Itreatment) 66min (with DNase Itreatment)







MD 757696



RBC Bioscience Corp. www.rbcbioscience.com info@rbcbioscience.com